

Tips for proper compression cap installation

1. Place the nut part of the compression cap on the tube and then slide the PTFE ferrule down shaft as shown in the picture below



2. Place the top part of the cap onto the tube so that the end of the tube rests as far inside the cap as possible as shown in the picture below.



3. Slide the nut along with the PTFE ferrule up the shaft and connect to the cap by screwing the 2 pieces together trapping the ferrule in between as shown in the picture below. **Hand tighten the compression cap and then use wrenches (9/16" & 1/2") to tighten another 1/4 turn. Gloves Recommended when handling tubes**

