



# OPERATIONAL INSTRUCTIONS FOR MILLIFLUIDIC DEVICE #FPTCMDO 9018\*

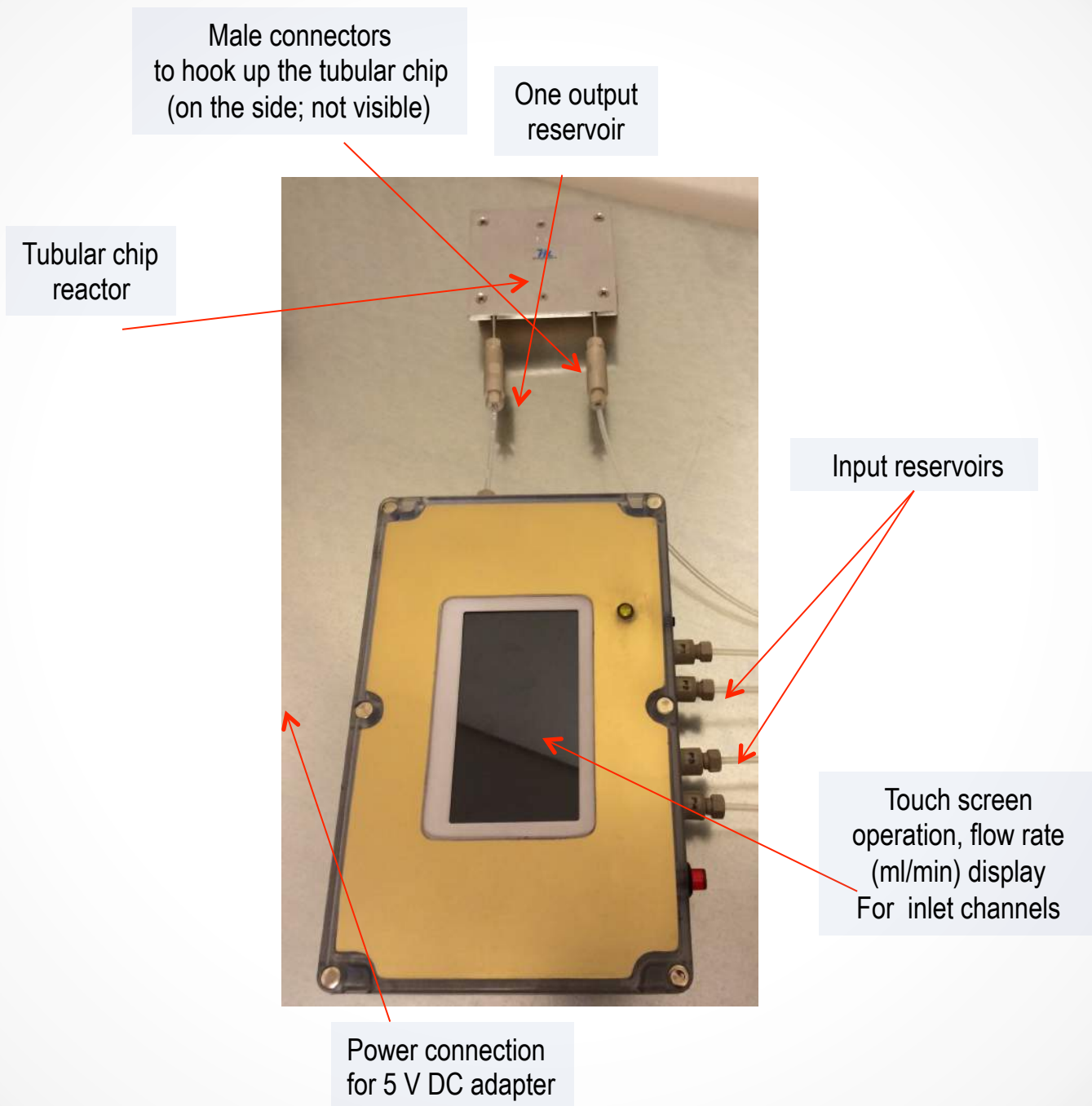


**\* DO NOT OPEN OR TAMPER WITH THE DEVICE. IN CASE OF  
ANY PROBLEM CONTACT THE CUSTOMER SERVICE  
IMMEDIATELY.**


# OPERATIONAL INSTRUCTIONS \*

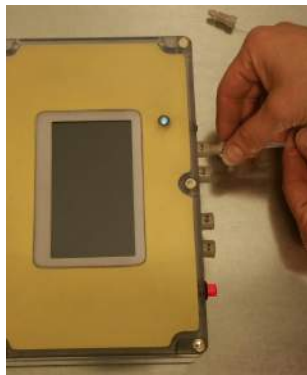


Step1a-connect  
input connectors 1-4

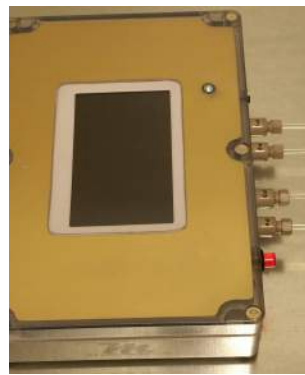


Step1b-1-4  
input  
connectors

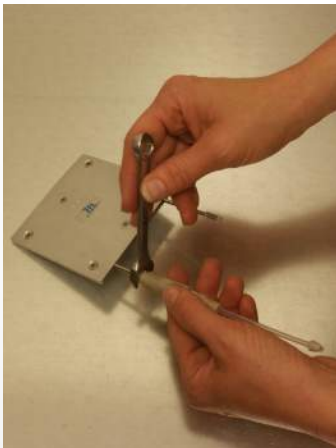

Step2-attach  
connectors to heat  
exchange

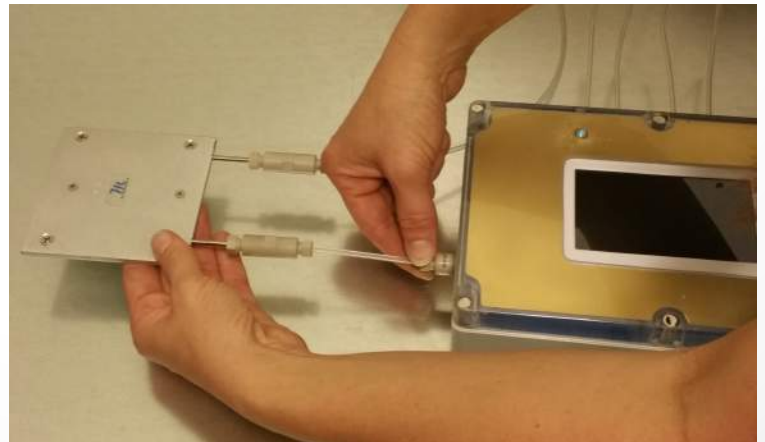

Step3-connect heat  
exchange to output

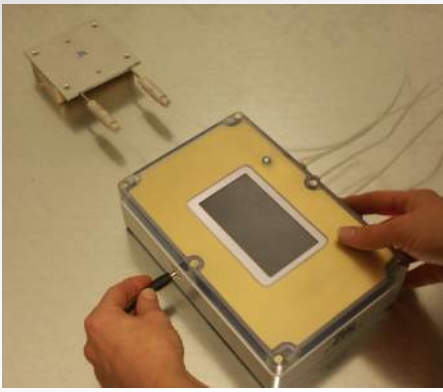


Step4-plug in power supply




Step5-turn on-touch screen




Step6-slide bar to right

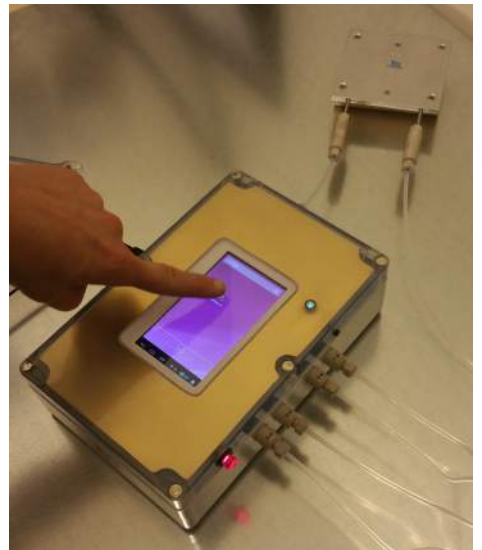

Step7-select Millifluidica app

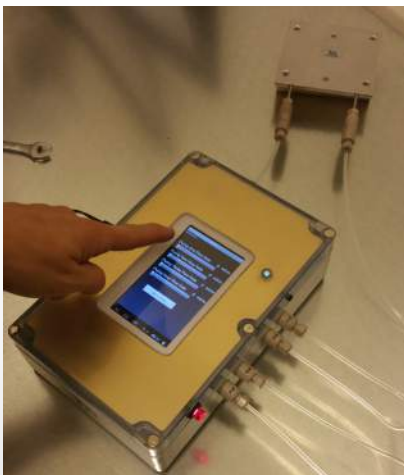


Step8-calibrate flow rates for pumps 1-4

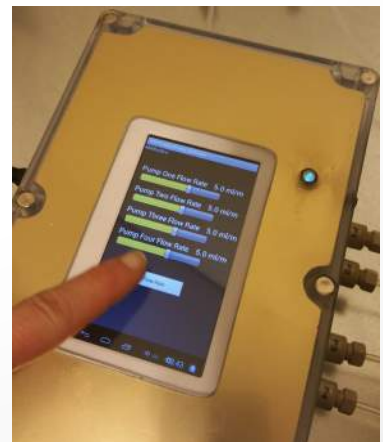

Step9-set flow rates



Step10-turn on pumps

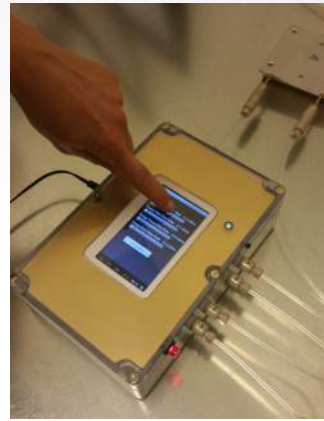

Step12-set flow rates

- The operator of the device is responsible for all actions and outcomes when using the device and agree to operate the device at his/her own risk. The manufacturer is not liable for any damages due to unanticipated accidents or misuse of the device or its components

**Questions? Call Customer Service: (650) 409 6546**